



CTR


SCIENTIFIC
www.ctr.com.mx

Versatile and Easy to Use



Thermo Scientific Nunc EasyFill Cell Factory

- Versatile
- Easy to use
- Bridges small scale research and commercial production
- Sterile, single use
- Reduces contamination risk

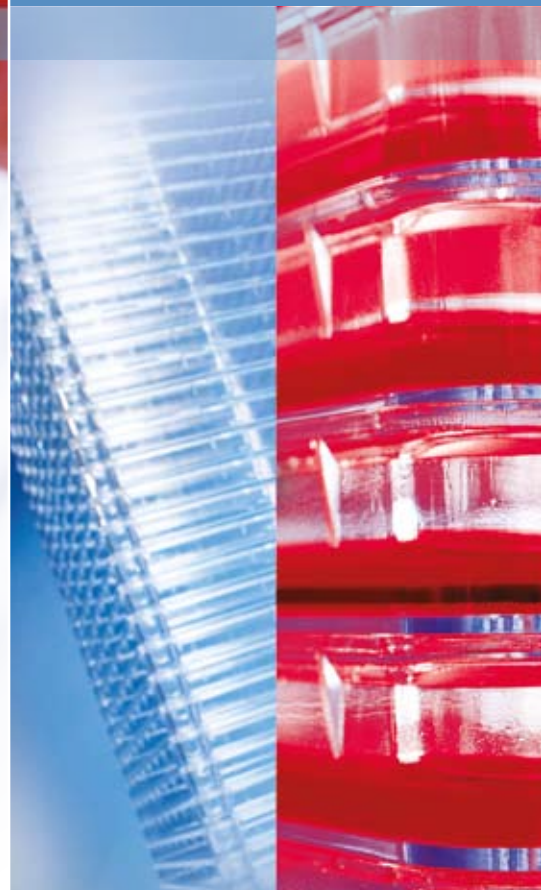

Instructions for use

Thermo Scientific Nunc EasyFill Cell Factory



1. Pour media directly into the Nunc EasyFill™ Cell Factory.



2. Equilibrate by placing the Cell Factory on the side with the small port.



3. Turn the Cell Factory 90°, so that the filling inlet is up. Medium will be separated with equal amounts in each chamber.



4. With filling inlet up, carefully tilt the Cell Factory to a horizontal incubation position. **Handle by holding on to the sides - do not manipulate by holding on to the upper rim.**



5. Incubate following appropriate protocol.



6. Loosen filter cap and remove cover cap. Empty directly into reservoir.



Tubing Configurator – Design your own integrated system

The online custom tubing set configurator lets you design your own integrated fluid transfer and bioprocess systems using Nalgene® and Nunc brand products. The easy-to-use browser interface enables users to choose from more than 150 standard components for fluid handling and generate

a two-dimensional CAD drawing to verify the design and assure its functionality.

For more information, go to www.tubingconfigurator.com

EasyFill Cell Factory Nunclon™ Δ

Polystyrene. Sterile
Length 335 mm, Width 205 mm



Cat. No.	140000	140250	140360	140400
Number of trays	1	2	4	10
Culture area, cm ²	630	1260	2520	6300
Suggested working vol., ml	200	400	800	2000
Units per pack/case	1/6	1/6	1/4	1/2

For scientific references visit our website at: www.nuncbrand.com/go/easyfill

Accessories

Cat. No.	173248	173249	140099	146008
Description	Gelman Filter with Connector	0.2 µm Air Filter with Connector	PC Connector	Filter Cap
Sterile	+	+	-	+
Units per pack/case	1/2	1/2	10/10	1/10

© 2008 Thermo Fisher Scientific Inc. All Rights Reserved. 76282/N20221 - Ver. 1.0 - 10/2008 - YNI.
All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.

North America - Tel: 1-800-625-4327 - technical.nunc@thermofisher.com
Asia Pacific - Tel: +65 6770 2807 - intlmtkg@thermofisher.com
China - Tel: 86-21-68654588 - info.nnchina@thermofisher.com
Europe - Tel: +45 4631 2000 - nunc.nalgene@thermofisher.com
India - Tel: +91-22-67162200 - Toll Free: 1800 22 8374 - contact.lpg.in@thermofisher.com
Japan - Tel: +81 3 3816 3355 - info@nalgenuc.co.jp - www.nalgenuc.co.jp
All other locations (USA, International Department) - Tel: +1 585 899 7198 - intlmtkg@thermofisher.com

www.thermo.com/easyfill

Thermo
SCIENTIFIC